

CARLOS ARBOLES S.A.

GRIFERÍA DE LABORATORIO / LABORATORY TAPS
 DUCHAS Y LAVAOJOS DE EMERGENCIA / EMERGENCY SHOWERS AND EYE WASH
 Compositor Narváez, 10 - Pol. Ind. Can Jardí - 08191 Rubí (Barcelona) - España
 Tel: +34 935 860 780 - Fax: +34 936 973 226
 info@carlosarboles.com - www.carlosarboles.com

TOLERANCIA GENERAL ISO 2768 - m
 TOLERANCE IN GENERAL ISO 2768 - m


ESCALA / SCALE
 1:20

DIBUJADO / DESIGNED: Xavier Mañes
 REVISADO / REVISED: Miguel Ángel V. H. 08/10/2010
 APROBADO / APPROVED: Xavier Mañes 08/10/2010
 REVISIÓN / REVISION: 1
 MATERIAL: -

MODELO / MODEL: 104220

DUCHA / LAVAOJOS DE EMERGENCIA PEDESTAL, RECOGEDOR / ROCIADOR ABS
 PEDESTAL EMERGENCY SHOWER / EYEWASH, ABS BOWL / SHOWER HEAD

*Este plano es propiedad de Carlos Arboles S.A. estando prohibida su reproducción y entrega a terceros sin previa autorización de Carlos Arboles S.A.
 This drawing is owned by Carlos Arboles S.A. prohibited its reproduction and delivery to third parties without prior permission of Carlos Arboles S.A.*

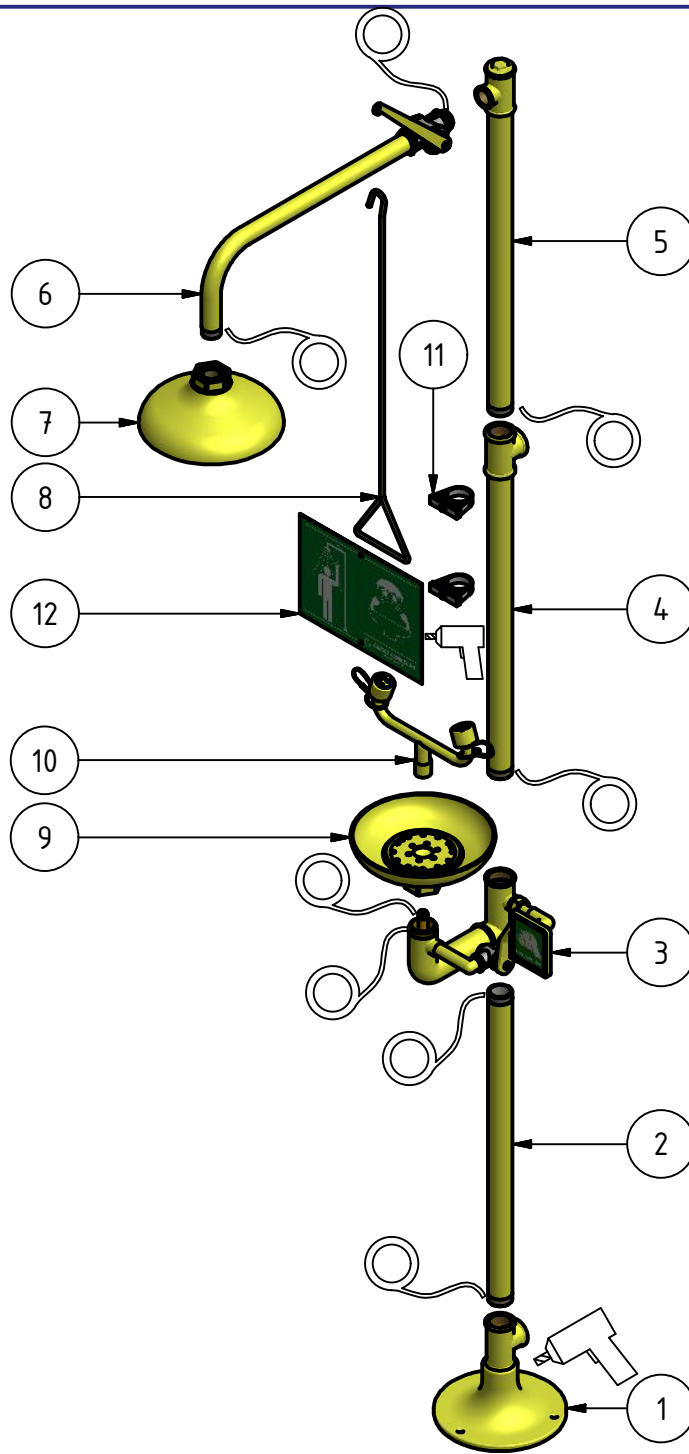

Lista de elementos / List of elements

Elemento / Element	Cantidad / Quantity	Modelo / Model	Descripción / Description
1	1	180030	CONJUNTO PIE / SUPPORT ASSEMBLY
2	1	864650	TUBO 1" 1/4 x 650 mm / PIPE 1" 1/4 x 650 mm
3	1	180135	CONJUNTO BLOQUE LAVAOJOS / EYEWASH BLOCK ASSEMBLY
4	1	180125	TUBO 1" 1/4 x 650 mm + TE 1" 1/4 / PIPE 1" 1/4 x 650 mm + T 1" 1/4
5	1	180120	TUBO 1" 1/4 x 650 mm + TE 1" 1/4 x 1" + TAPÓN / PIPE 1" 1/4 x 650 mm + T 1" 1/4 x 1" + PLUG
6	1	180075	CONJUNTO BRAZO DUCHA / SHOWER ARM ASSEMBLY
7	1	190100	ROCIADOR DUCHA ABS / ABS SHOWER HEAD
8	1	190105	TIRADOR DUCHA / PULL ROD
9	1	190150	RECOGEDOR LAVAOJOS ABS / ABS BOWL
10	1	190136	DISTRIBUIDOR LAVAOJOS / EYEWASH ASSEMBLY
11	1	191007	KIT FIJACIÓN PARA LETRERO / FIXING KIT
12	1	191006	LETRERO DUCHA/LAVAOJOS / SHOWER/EYEWASH SIGN



CARLOS ARBOLES S.A.

GRIFERÍA DE LABORATORIO / LABORATORY TAPS
 DUCHAS Y LAVAOJOS DE EMERGENCIA / EMERGENCY SHOWERS AND EYE WASH
 Compositor Narváez, 10 - Pol. Ind. Can Jardí - 08191 Rubí (Barcelona) - España
 Tel: +34 935 860 780 - Fax: +34 936 973 226
 info@carlosarboles.com - www.carlosarboles.com

TOLERANCIA GENERAL ISO 2768 - m
 TOLERANCE IN GENERAL ISO 2768 - m

ESCALA / SCALE
 1:13

DIBUJADO / DESIGNED: Xavier Mañes
 REVISADO / REVISED: Miguel Ángel V. H. 08/10/2010
 APROBADO / APPROVED: Xavier Mañes 08/10/2010
 REVISIÓN / REVISION: 1
 MATERIAL: -

MODELO / MODEL: 104220

DUCHA / LAVAOJOS DE EMERGENCIA PEDESTAL, RECOGEDOR / ROCIADOR ABS
 PEDESTAL EMERGENCY SHOWER / EYEWASH, ABS BOWL / SHOWER HEAD